

Instructions


The touch pad is the whole grey numpad area. Toggle off to use 'CTRL-ALT-DEL' to login, toggle on again to use as a mouse.

- Left click with a one finger tap.
- Right click with a two finger tap.
- Left click drag (to move a window, select items or do phase-correction in Topspin) use three fingers and drag.
- Scroll with a two finger drag.

Once you have finished using it, hold the lock key (3) for 3 seconds to disable the keyboard and wipe clean with the sanitising fluid.

These keyboards are good for simple operations, complicated processing should be done on your own computer.

Remember that CTRL-Z will undo the last operation!

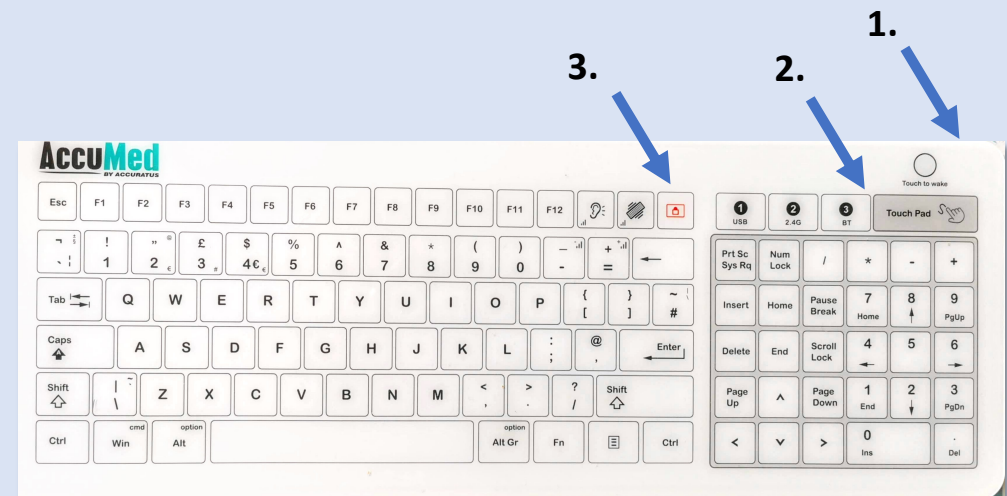

1 – Power button is on the back.

2 – Toggles touch pad on/off

3 – Keyboard lock. Hold to engage/release lock.

Any problems, please email nmr@ch.cam.ac.uk